


Revolutionize Your Research

Leading researchers are discovering the promise of the SELECT SERIES™ Cyclic™ IMS. This revolutionary new instrument combines novel cyclic ion mobility separation with state-of-the-art, high-performance, time-of-flight mass spectrometry (ToF), to help bring you to the forefront of scientific discovery.

Discover a New Path

The circular path of the SELECT SERIES Cyclic

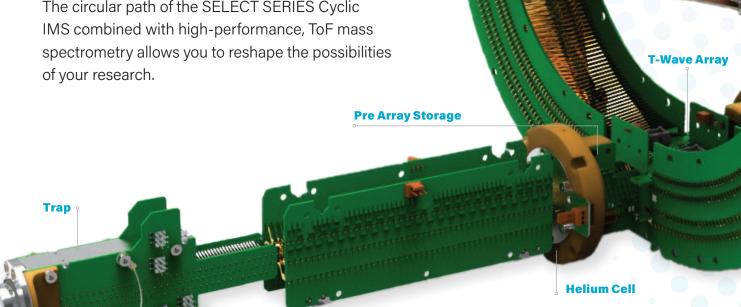

Combines cyclic ion mobility (cIM) separation with ToF mass spectrometry

Separates more than ever before with ultimate ion mobility resolution for complete confidence

Performs unique IMSⁿ experiments for increased selectivity and advanced characterization

Enables premobility mass selection and mobility selection of precursor and product ions


Provides three distinct fragmentation regions, enabling increased experimental flexibility for solving complex problems

Cyclic IMS Ion Optics





Now you can acquire a wide range of experiments, including single and multi-pass ion mobility, IMS/IMS, and IMSⁿ, all combined with high resolution ToF data acquisition.

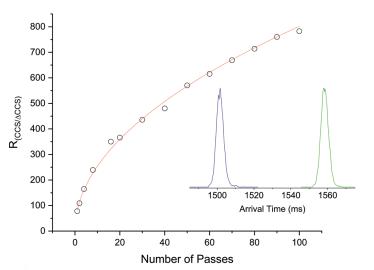

XS Transfer Device

ULTIMATE MOBILITY RESOLUTION

IMS resolution increases with the number of passes around the cIM device.
Demonstrated here by the separation of reverse sequence peptides SDGRG and GRGDS. The Arrival Time Distributions (ATDs) obtained after 100 passes around the device are inset.

NOVEL IMSⁿ EXPERIMENTS

The unique instrument geometry and software design enables you to perform a wide range of experiments such as: ion selection of specific mobility separated species, ion storage, activation, multiple rounds of IMS/IMS (IMSⁿ), and importantly, custom combinations. Experiments can be designed and implemented with varying storage and separation times to provide bespoke methods to investigate structural and positional isomers, gas phase interconversion of molecular structures, and ion conformations.


SUPERIOR TOF PERFORMANCE

Have complete confidence in your results with improved sensitivity, resolution >100,000 (FWHM), reliable exact mass measurements (low ppm), and improved dynamic range.

REMARKABLE TOF RESOLUTION

Mass spectrum highlighting resolution of >100,000 (FWHM) on the 7+ charge state of monomeric streptavidin.



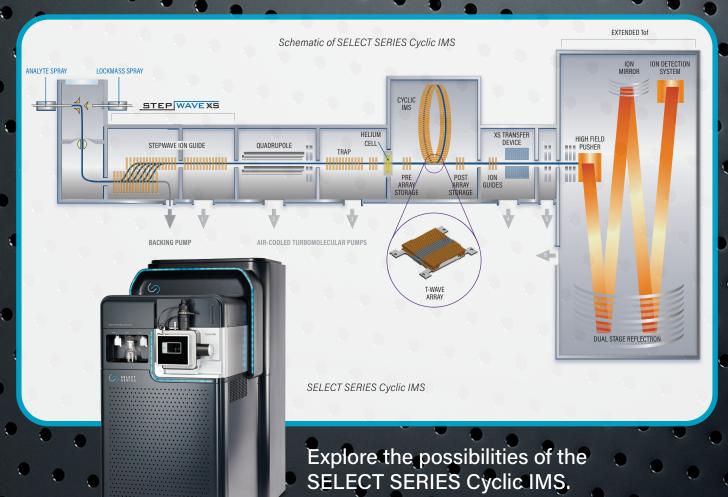

For your local sales office, please visit waters.com/contact



Waters Corporation 34 Maple Street Milford, MA 01757 U.S.A. T: 1 508 478 2000 F: 1 508 872 1990 www.waters.com

Waters

www.waters.com/CyclicIMS

THE SCIENCE OF WHAT'S POSSIBLE."

Waters, The Science of What's Possible, SELECT SERIES, and Cyclic are trademarks of Waters Corporation. All other trademarks are the property of their respective owners.

©2020 Waters Corporation. Produced in the U.S.A. March 2020 720006752EN TC-SIG